

## Setup

Place Pipette Shaker on a flat, stable surface. Ensure that the controls are set in the “off” position. Plug into the appropriate electrical outlet.

## Controls

Three-position Power Switch: Used to turn the shaker on and off. Positions are “ON,” “Off,” and “Timer.” “On” setting runs continuously; “Timer” setting runs for periods set on timer.

Red Power Light: Signals that unit is on

## Operation


- 1 Place samples in the top holder.
- 2 Flip the Power Switch to either the “On” or “Timer” position.

## Specifications

Specifications:  
Speed: 2500 RPM  
110V / 60Hz (220V available)  
Height: 158.75mm (6.25 in)  
Width: 133.35mm (5.25 in)  
Weight: 1.9Kg (4.2 lbs.)



Notes for Safe Use:  
Do not operate unit in or near water or other liquids.  
Unplug unit before cleaning or repairing.  
Do not use if electrical cord is damaged.  
Replace cord if damaged.

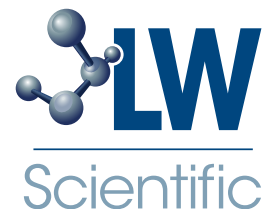

# Pipette Shaker

Instruction Manual



Models pictured:  
Pipette Shaker (SHL-PPF7-06F1)

See back page for model specifications.



## Unpacking

Carefully unpack the pipette shaker, saving packing materials in the unlikely event you need to return the rocker. Immediately call LW Scientific if you detect any damage. If you need to return the turbo mixer, please use the original packing material.