


Turbo Mixer

Instruction Manual

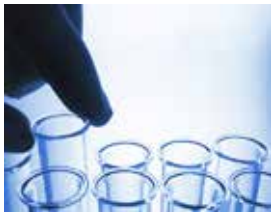

Models pictured:
MXL-TMV7-7RC1

See back page for model specifications.



Unpacking

Carefully unpack the Turbo Mixer, saving packing materials in the unlikely event you need to return the mixer. Immediately call LW Scientific if you detect any damage. If you need to return the turbo mixer, please use the original packing material.



Note: Always treat specimen samples as though they are infectious. Do not autoclave.

Setup

Place Turbo Mixer on a flat, stable surface. Ensure that the controls are set in the "off" position. Plug into the appropriate electrical outlet.

Controls

Three-position Power Switch: Used to turn the blood rocker on and off. Positions are "Auto," "Off," and "Touch." "Auto" setting runs continuously; "Touch" setting runs only when items are placed on the platform or in the cup.

Speed Control Knob: Used to control variable speed, up to 3000 rpm

Operation

- 1 Place samples to be mixed on the top of the platform or in the cup.
- 2 Flip the Power Switch to either the "Auto" or "Touch" position.

Specifications

Speed:	3000 RPM
Power:	110V / 60Hz (220V available)
Width:	160mm (6.3")
Depth:	130mm (5.2")
Height:	170mm (6.7")
Weight:	3.2Kg (7lbs)



Notes for Safe Use:

- Do not operate unit in or near water or other liquids.
- Unplug unit before cleaning or repairing.